APPENDIX J

RECOMMENDED PROCEDURES FOR PREPARING WHOLE FISH COMPOSITE HOMOGENATE SAMPLES

APPENDIX J

RECOMMENDED PROCEDURES FOR PREPARING WHOLE FISH COMPOSITE HOMOGENATE SAMPLES

J.1 GENERAL GUIDELINES


Laboratory processing to prepare whole fish composite samples (diagrammed in Figure J-1) involves

- Inspecting individual fish for foreign material on the surface and rinsing if necessary
- Weighing individual fish
- Examining each fish for morphological abnormalities (optional)
- Removing scales or otoliths for age determination (optional)
- Determining the sex of each fish (optional)
- Preparing individual whole fish homogenates
- Preparing a composite whole fish homogenate.


Whole fish should be shipped on wet or blue ice from the field to the sample processing laboratory if next-day delivery is assured. Fish samples arriving in this manner (chilled but not frozen) should be weighed, scales and/or otoliths removed, and the sex of each fish determined within 48 hours of sample collection. The grinding/homogenization procedure may be carried out more easily and efficiently if the sample has been frozen previously (Stober, 1991). Therefore, the samples should then be frozen (\leq -20 °C) in the laboratory prior to being homogenized.

If the fish samples arrive frozen (i.e., on dry ice) at the sample processing laboratory, precautions should be taken during weighing, removal of scales and/or otoliths, and sex determination to ensure that any liquid formed in thawing remains with the sample. **Note:** The liquid will contain target analyte contaminants and lipid material that should be included in the sample for analysis.

The thawed or partially thawed whole fish should then be homogenized individually, and equal weights of each homogenate should be combined to form the composite sample. Individual homogenates and/or composite homogenates may be frozen; however, frozen individual homogenates must be rehomogenized before compositing, and frozen composite homogenates must be rehomogenized before aliquotting for analysis. The maximum holding time from sample collection to analysis for mercury is 28 days at \leq -20 °C; for all other analytes, the holding time is 1 year at \leq -20 °C (Stober, 1991). Recommended container materials,

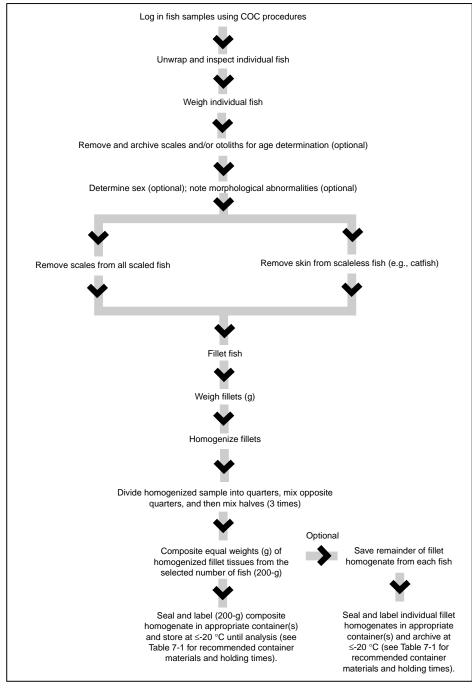

COC = Chain of custody.

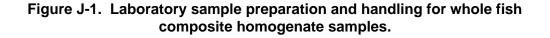

preservation temperatures, and holding times are given in Table J-1. **Note:** Holding times in Table J-1 are maximum times recommended for holding samples from the time they are received at the laboratory until they are analyzed. These holding times are based on guidance that is sometimes administrative rather than technical in nature; there are no promulgated holding time criteria for tissues (U.S. EPA, 1995b). If states choose to use longer holding times, they must demonstrate and document the stability of the target analyte residues over the extended holding times.

J.2 SAMPLE PROCESSING PROCEDURES

Fish sample processing procedures are discussed in more detail in the sections below. Each time custody of a sample or set of samples is transferred from one person to another during processing, the Personal Custody Record of the chainof-custody (COC) form that originated in the field (Figure 6-8) must be completed and signed by both parties so that possession and location of the samples can be traced at all times (see Section 7.1). As each sample processing procedure is performed, it should be documented directly in a bound laboratory notebook or on standard forms that can be taped or pasted into the notebook. The use of a standard form is recommended to ensure consistency and completeness of the record. Several existing programs have developed forms similar to the sample processing record for whole fish composite samples shown in Figure J-2.

J.2.1 Sample Inspection

Individual fish received for filleting should be unwrapped and inspected carefully to ensure that they have not been compromised in any way (i.e., not properly preserved during shipment). Any specimen deemed unsuitable for further processing and analysis should be discarded and identified on the sample processing record.

J.2.2 Sample Weighing

A wet weight should be determined for each fish. All samples should be weighed on balances that are properly calibrated and of adequate accuracy and precision to meet program data quality objectives. Balance calibration should be checked at the beginning and end of each weighing session and after every 20 weighings in a weighing session.

Fish shipped on wet or blue ice should be weighed directly on a foil-lined balance tray. To prevent cross contamination between individual fish, the foil lining should be replaced after each weighing. Frozen fish (i.e., those shipped on dry ice) should be weighed in clean, tared, noncontaminating containers if they will thaw before the weighing can be completed. Liquid from the thawed sample must be

	Matrix	Sample container	Storage		
Analyte			Preservation	Holding time ^a	
Mercury	Tissue (whole specimens, homogenates)	Plastic, borosilicate glass, quartz, and PTFE	Freeze at <u><</u> -20 °C	28 days⁵	
Other metals	Tissue (whole specimens, homogenates)	Plastic, borosilicate glass, quartz, and PTFE	Freeze at <u><</u> -20 °C	6 months ^c	
Organics	Tissue (whole specimens, homogenates)	Borosilicate glass, quartz, PTFE, and aluminum foil	Freeze at <u><</u> -20 °C	1 year ^d	
Metals and organics	Tissue (whole specimens, homogenates)	Borosilicate glass, quartz, and PTFE	Freeze at <u><</u> -20 °C	28 days (mercury); 6 months (for other metals); and 1 year (for organics)	
Lipids	Tissue (whole specimens, homogenates)	Plastic, borosilicate glass, quartz, PTFE	Freeze at <u><</u> -20 °C	1 year	

Table J-1. Recommendations for Container Materials, Preservation, and HoldingTimes for Fish, Shellfish, and Turtle Tissues from Receipt at Sample ProcessingLaboratory to Analysis

PTFE = Polytetrafluoroethylene for Teflon.


^a Maximum holding times recommended by U.S. EPA (1995b).

^d This maximum holding time is also recommended by the Puget Sound Estuary Program (1990). The California Department of Fish and Game (1990) and the USGS National Water Quality Assessment Program (Crawford and Luoma, 1993) recommend a more conservative maximum holding time of 6 months. EPA (1995a) recommends a maximum holding time of 1 year at ≤-10 °C for dioxins and dibenzofurans.

^b This maximum holding time is also recommended by the Puget Sound Estuary Program (1990). The California Department of Fish and Game (1990) and the USGS National Water Quality Assessment Program (Crawford and Luoma, 1993) recommend a maximum holding time of 6 months for all metals, including mercury.

^c This maximum holding time is also recommended by the California Department of Fish and Game (1990), the 301(h) monitoring program (U.S. EPA, 1986), and the USGS National Water Quality Assessment Program (Crawford and Luoma, 1993). The Puget Sound Estuary Program (1990) recommends a maximum holding time of 2 years.

Project No.			_ Sampling Date and Time:				
SITE LOCA		g 🗋 ;	Intens	ive: Phase I	Phase II		
				Lat / ong ·			
•		umber:					
	-	lame:					
Composite Sample #:							
Fish # 001	Weight (g)	Scales/Otoliths Removed (-⁄)	Sex (M, F)	Homogenate Prepared (✔)	Weight of homogenate taken for composite (g		
002 _							
003 _					<u></u>		
004 _	-						
005 _		- <u></u>					
006 _							
007 _							
008 _							
009 _							
010 _ Analyst							
Initials/Date _	/	/	/	/	/		
			Total Com	posite Homogenate \	Weight		
Predator –	Species Name:				· · · · · · · · · · · · · · · · · · ·		
Composite	Sample #:		Number of Individuals:				
Fish #	Weight (g)	Scales/Otoliths Removed (✓)	Sex (M, F)	Homogenate Prepared (✔)	Weight of homogenate taken for composite (g		
001 002							
002							
004			—	and an and a second			
005							
006 _							
007 _							
008 _							
009 _		<u></u>					
010 _							
Analyst Initials/Date _	/	/	/	/	/		
		Total Composite Homogenate Weight					

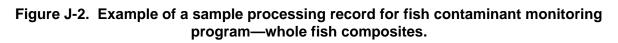

kept in the container as part of the sample because it will contain lipid material that has separated from the tissue (Stober, 1991).

All weights should be recorded to the nearest gram on the sample processing record and/or in the laboratory notebook.

J.2.3 Age Determination

Age provides a good indication of the duration of exposure to pollutants (Versar, 1982). A few scales or otoliths (Jearld, 1983) should be removed from each fish and delivered to a fisheries biologist for age determination. For most warm water inland gamefish, 5 to 10 scales should be removed from below the lateral line and behind the pectoral fin. On softrayed fish such as trout and salmon, the scales should be taken just above the lateral line (WDNR, 1988). For catfish and other scaleless fish, the pectoral fin spines should be clipped and saved (Versar, 1982). The scales, spines, or otoliths may be stored by sealing them in small envelopes (such as coin envelopes) or plastic bags labeled with, and cross-referenced by, the identification number assigned to the tissue specimen (Versar, 1982). Removal of scales, spines, or otoliths from each fish should be noted (by a check mark) on the sample processing record.

J.2.4 Sex Determination (Optional)

To determine the sex of a fish, an incision should be made on the ventral surface of the body from a point immediately anterior to the anus toward the head to a point immediately posterior to the pelvic fins. If necessary, a second incision should be made on the left side of the fish from the initial point of the first incision toward the dorsal fin. The resulting flap should be folded back to observe the gonads. Ovaries appear whitish to greenish to golden brown and have a granular texture. Testes appear creamy white and have a smooth texture (Texas Water Commission, 1990). The sex of each fish should be recorded on the sample processing record.

J.2.5 Assessment of Morphological Abnormalities (Optional)

Assessment of gross morphological abnormalities in finfish is optional. This assessment may be conducted in the field (see Section 6.3.1.5) or during initial inspection at the central processing laboratory prior to filleting. States interested in documenting morphological abnormalities should consult Sinderman (1983) and review recommended protocols for fish pathology studies used in the Puget Sound Estuary Program (1990).

J.2.6 Preparation of Individual Homogenates

To ensure even distribution of contaminants throughout tissue samples, whole fish must be ground and homogenized prior to analyses.

Smaller whole fish may be ground in a hand crank meat grinder (fish < 300 g) or a food processor (fish 300-1,000 g). Larger (>1,000 g) fish may be cut into 2.5-cm cubes with a food service band saw and then ground in either a small or large homogenizer. To avoid contamination by metals, grinders and homogenizers used to grind and blend tissue should have tantalum or titanium blades and/or probes. Stainless steel blades and probes have been found to be a potential source of nickel and chromium contamination (due to abrasion at high speeds) and should be avoided.

Grinding and homogenization of biological tissue, especially skin from whole fish samples, is easier when the tissue is partially frozen (Stober, 1991). Chilling the grinder/homogenizer briefly with a few chips of dry ice will reduce the tendency of the tissue to stick to the grinder.

The ground sample should be divided into quarters, opposite quarters mixed together by hand, and the two halves mixed back together. The grinding, quartering, and hand mixing should be repeated two more times. If chunks of tissue are present at this point, the grinding/homogenizing should be repeated. No chunks of tissue should remain because these may not be extracted or digested efficiently. If the sample is to be analyzed for metals only, the ground tissue may be mixed by hand in a polyethylene bag (Stober, 1991). Homogenization of each individual fish should be noted on the sample processing record. At this time, individual whole fish homogenates may be either composited or frozen and stored at \leq -20 °C in cleaned containers that are noncontaminating for the analyses to be performed (see Table J-1).

J.2.7 Preparation of Composite Homogenates

Composite homogenates should be prepared from equal weights of individual homogenates. If individual whole fish homogenates have been frozen, they should be thawed partially and rehomogenized prior to compositing. Any associated liquid should be maintained as a part of the sample. The weight of each individual homogenate that is used in the composite homogenate should be recorded, to the nearest gram, on the sample processing record.

Each composite homogenate should be blended by dividing it into quarters, mixing opposite quarters together by hand, and mixing the two halves together. The quartering and mixing should be repeated at least two more times. If the sample is to be analyzed only for metals, the composite homogenate may be mixed by hand in a polyethylene bag (Stober, 1991). At this time, the composite homogenate may be processed for analysis or frozen and stored at \leq -20 °C (see Table J-1).

The remainder of each individual homogenate should be archived at \leq -20 °C with the designation "Archive" and the expiration date recorded on the sample label. The location of the archived samples should be indicated on the sample processing record under "Notes."

It is essential that the weights of individual homogenates yield a composite homogenate of adequate size to perform all necessary analyses. Weights of individual homogenates required for a composite homogenate, based on the number of fish per composite and the weight of composite homogenate recommended for analyses of all screening study target analytes (see Table 4-1), are given in Table J-2. The total composite weight required for intensive studies may be less than in screening studies if the number of target analytes is reduced significantly.

The recommended sample size of 200 g for screening studies is intended to provide sufficient sample material to (1) analyze for all recommended target analytes (see Table 4-1) at appropriate detection limits, (2) meet minimum QA and QC requirements for the analyses of replicate, matrix spike, and duplicate matrix spike samples (see Section 8.3.3.4), and (3) allow for reanalysis if the QA and QC control limits are not met or if the sample is lost. However, sample size requirements may vary among laboratories and the analytical methods used. Therefore, it is the responsibility of each program manager to consult with the analytical laboratory supervisor to determine the actual weights of composite homogenates required to analyze for all selected target analytes at appropriate detection limits.

J.3 REFERENCES

- California Department of Fish and Game. 1990. *Laboratory Quality Assurance Program Plan.* Environmental Services Division, Sacramento, CA.
- Crawford, J.K., and S.N. Luoma. 1993. *Guidelines for Studies of Contaminants in Biological Tissues for the National Water-Quality Assessment Program.* USGS Open-File Report 92-494. U.S. Geological Survey, Lemoyne, PA.
- Jearld, A. 1983. Age determination. pp. 301-324. In: *Fisheries Techniques*. L.A. Nielsen and D. Johnson (eds.). American Fisheries Society, Bethesda, MD.
- Puget Sound Estuary Program. 1990 (revised). Recommended protocols for fish pathology studies in Puget Sound. Prepared by PTI Environmental Services, Bellevue, WA. In: *Recommended Protocols and Guidelines for Measuring Selected Environmental Variables in Puget Sound.* Region 10, U.S. Environmental Protection Agency, Seattle, WA. (Looseleaf)
- Sinderman, C. J. 1983. An examination of some relationships between pollution and disease. *Rapp. P. V. Reun. Cons. Int. Explor. Mer.* 182:37-43.
- Stober, Q. J. 1991. Guidelines for Fish Sampling and Tissue Preparation for Bioaccumulative Contaminants. Environmental Services Division, Region 4, U.S. Environmental Protection Agency, Athens, GA.

- Texas Water Commission. 1990. *Texas Tissue Sampling Guidelines.* Texas Water Commission, Austin, TX.
- U.S. EPA (U.S. Environmental Protection Agency). 1986. Bioaccumulation Monitoring Guidance: 4. Analytical Methods for U.S. EPA Priority Pollutants and 301(h) Pesticides in Tissues from Marine and Estuarine Organisms. EPA-503/6-90-002. Office of Marine and Estuarine Protection, Washington, DC.
- U.S. EPA (U.S. Environmental Protection Agency). 1995a. *Method 1613b. Tetra- through Octa-Chlorinated Dioxins and Furans by Isotope Dilution HRGC/HRMS.* Final Draft. Office of Water, Office of Science and Technology, Washington, DC.
- U.S. EPA (Environmental Protection Agency). 1995b. QA/QC Guidance for Sampling and Analysis of Sediments, Water, and Tissues for Dredged Material Evaluations—Chemical Evaluations. EPA 823-B-95-001. Office of Water, Washington, DC, and Department of the Army, U.S. Army Corps of Engineers, Washington, DC.
- Versar, Inc. 1982. Sampling Protocols for Collecting Surface Water, Bed Sediment, Bivalves and Fish for Priority Pollutant Analysis--Final Draft Report. EPA Contract 68-01-6195. Prepared for U.S. EPA Office of Water Regulations and Standards. Versar, Inc., Springfield, VA.
- WDNR (Wisconsin Department of Natural Resources). 1988. Fish Contaminant Monitoring Program—Field and Laboratory Guidelines (1005.1). Madison, WI.